

*P. heparinus* Culture Medium recipes

1. 1% yeast broth (1 g yeast extract/100 ml dH<sub>2</sub>O. Autoclave.)
2. Non-spreading plates (1.0% yeast extract and 1.5% agar; For 100 ml: 1 g yeast extract, 1.5 g agar and water to 100 ml. Autoclave.)
3. Spreading plates (0.5% yeast extract and .5% agarose; For 100 ml: 0.5 g yeast extract, .05 g agarose and water to 100 ml. Autoclave.)